

user manual

# **pco.**firmware upgrader pco.edge USB 3.0

1288   
EMVA Standard Compliant

**pco.**  
An Excelitas Technologies Brand

**PCO asks you to carefully read and follow the instructions in this document.  
For any questions or comments, please feel free to contact us at any time.**




An Excelitas Technologies Brand

telephone:	+49 (0) 9441 2005 50
fax:	+49 (0) 9441 2005 20
postal address:	Excelitas PCO GmbH Donaupark 11 93309 Kelheim, Germany
email:	<a href="mailto:info@pco.de">info@pco.de</a>
web:	<a href="http://www.pco.de">www.pco.de</a>

pco.firmware upgrader pco.edge USB 3.0 user manual 1.5.0

Released March 2022

©Copyright Excelitas PCO GmbH

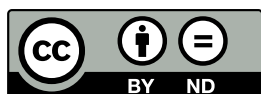

This work is licensed under the Creative Commons Attribution-NoDerivatives 4.0 International License. To view a copy of this license, visit <http://creativecommons.org/licenses/by-nd/4.0/> or send a letter to Creative Commons, PO Box 1866, Mountain View, CA 94042, USA.

# Contents

<b>1</b>	<b>General</b>	<b>4</b>
1.1	Introduction . . . . .	4
1.2	Availability . . . . .	4
1.3	Support . . . . .	4
<b>2</b>	<b>Firmware upgrade preparations</b>	<b>5</b>
<b>3</b>	<b>Upgrade process</b>	<b>6</b>

# 1 General

## 1.1 Introduction

This manual describes the firmware upgrade procedure of specific PCO cameras (see **1.2**).

There are different versions of the firmware upgrader for 32 bit and 64 bit computer systems.

## 1.2 Availability

Available for these camera models:

- pco.edge 5.5, pco.edge 4.2, pco.edge 3.1 with USB 3.0 interface

**Note** For all firmware versions from 1.06 to 1.23 and version 2.02.

## 1.3 Support

In case of errors please contact us at PCO:

- [support@pco.de](mailto:support@pco.de)
- tel +49 9441 2005 0



## 2 Firmware upgrade preparations

The following files are required for the firmware upgrade process:

- Firmware upgrader application: programmer\_pco.edge\_usb3\_upgrade.exe
- Firmware upgrade file (example): FW\_pco.edge\_4.2\_usb3.0\_203\_UPGRADE.ppg

Follow the steps below:

- 1 The camera must be connected to your computer.**
- 2 Now switch on the camera.**
- 3 Please close all open programs on your computer.**
- 4 Start the programmer\_pco.edge\_usb3\_upgrade.exe**

**Note** Do not unplug and/or switch off any of your devices, neither the computer, nor the camera during the update!

## 3 Upgrade process

Press **Load Firmware** and select the appropriate upgrade file.



If a wrong camera is connected or a wrong package is selected, the firmware upgrader will show a warning message.




If the firmware package and camera fit: The message **Firmware packet file is loaded successfully** will appear, then press **Program Camera**.



The upgrade process will take a few minutes. The camera system will restart during this process and the connection will be interrupted for a few minutes.



Wait until the message **PCO Camera Firmware Update succeeded. Close Application?** appears.

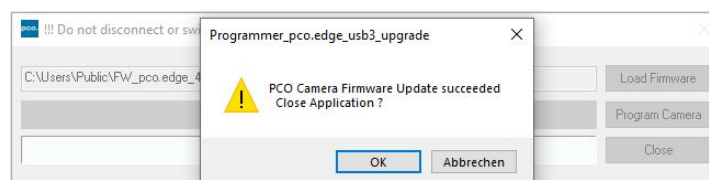

The firmware upgrade process is now finished. Press **OK**.

## contact

### pco europe

+49 9441 2005 50  
info@pco.de  
pco.de

### pco america

+1 866 678 4566  
info@pco-tech.com  
pco-tech.com

### pco asia

+65 6549 7054  
info@pco-imaging.com  
pco-imaging.com

### pco china

+86 512 67634643  
info@pco.cn  
pco.cn



for application stories  
please visit our website

# pco.

An Excelitas Technologies Brand